

BR100-C1-7L Microbial Fermentation System

Technical Specifications

Prepared by Labfirst Scientific, October 2023
 Email: sales@lab1st.com; Tel: +86-13524020331



1. General Information

Items Included	Control Console	BR100-C1
	Culture Vessel	7L Boro3.3 glass vessel with top mechanical stirring (25~75% working volume), SS316L cover
	Sensors	PT100, DO, foam, pH
	Feeding Pumps	4 peristaltic pump heads, controlled
	Gas Inlet	<ul style="list-style-type: none"> · Sparger from the bottom (Air + O₂ + N₂ + CO₂): controlled by rotameter and electromagnetic valve; · Overlay (Air): manually controlled by rotameter;
	Heating & Cooling	Heated by bottom electric hot plate; cooled by stainless steel cooling tube (for chilled water)
	Accessories	4 supplement bottles, SS316L feeding needle, etc.
	Consumables	O-rings, fuses, silicone stoppers, inlet air filter, vent filter, 20m silicone tubing, etc.
	Sensors Included	pH
DO		Sterilizable probe by Hamilton / Mettler Brand, with 1 meter electrode wire
Temperature		German made PT100 sensor, sterilizable
Foam		Foam probe, made in China

Trust us with your projects!

Lab1st has produced and sold more than 11,000 bioreactors, including 4,500+ Benchtop bioreactors and 6,000+ stainless steel fermenters since 1998. [3 production centers, 20+ years experience]

2. Specific Information

Functions Included	Temperature	<ul style="list-style-type: none"> · Robust PID algorithm; · Electrical-heating by SS316L electric hot plate at vessel bottom, chilled by tap or chilled water supply; · Sequential control coupled with fermentation process;
	pH	<ul style="list-style-type: none"> · pH control coupled with peristaltic pump or CO₂; · Manual pH control by peristaltic pump;
	DO	<ul style="list-style-type: none"> · Cascade control, coupled with different parameters (gas inlet and peristaltic pump);
	Agitation	<ul style="list-style-type: none"> · Robust PID algorithm; · Manual speed control or speed control coupled with other parameters;
	Foam	<ul style="list-style-type: none"> · Foam control coupled with peristaltic pump; (anti-foam); · Mechanical defoaming blade;
	Feeding (peristaltic pump)	<ul style="list-style-type: none"> · Accumulation volume control; · Manual control; Or control by plan or pre-set curve; · Coupled with other parameters (DO, pH, etc);
	Fermentation Sequence	Pre-set fermentation sequences along with timeline (pH, DO, agitation, temperature, etc)
	Others	<ul style="list-style-type: none"> · Historical data export and curves; · Fermentation/culture batch record; · Calibration of sensors and peristaltic pumps; · 3-level authority setting & control; · Record historical operations

2.1 Control Console

General Info	Model	BR100-C1
	Description	Desktop control console for cell culture fermentation, 1 console for 1 vessel
	Dimension	380×500×730mm
	Net Weight	60 Kg
	Touch Screen	PLC based 10" touch screen
	Communication	USB (data export), , Ethernet interface, RS485
	Electricity	208V-240V, 50-60Hz, single phase, 1.8 kW
Gas Inlet	Control	<ul style="list-style-type: none"> · Sparger from the bottom (Air + O₂ + N₂ + CO₂): controlled by rotameter and electromagnetic valve; · Overlay (Air): manually controlled by rotameter;
	Flow Rate	2 vvm
	Filtration	0.2 um
	Connection	8mm hose barb
Temperature	Range	8C above coolant to 40C above ambient (0-65C absolute)
	Control Accuracy	+/- 0.1C

Trust us with your projects!

Lab1st has produced and sold more than 11,000 bioreactors, including 4,500+ Benchtop bioreactors and 6,000+ stainless steel fermenters since 1998. [3 production centers, 20+ years experience]

pH	Range	2.00~12.00 (resolution 0.01)
	Control Accuracy	+/- 0.1
DO	Range	0~150% (resolution 0.1%)
	Control Accuracy	+/- 3%
Upgrades	Analysis or Control	OUR, CER, KLa, RQ,ORP, methanol, exhaust gas (O2, CO2), glucose, ect.
	Gas Supply Control	TMFC control (flow rate - automatic)
	Peristaltic Pump	More peristaltic pumps for different functions
	Scale	More peristaltic pumps for different functions; Upgrade to Watson-Marlow peristaltic pump heads;
	Light	Light source can be selected from red, blue, and white. Its intensity is adjustable [0-100%]

2.2 Vessel Information

General Info	Volume	7L max, with 25-75% working volume
	Material	Boro3.3 glass and SS316L (polishing better than Ra 0.4 um)
	Ratio	Diameter to height, about 1:2
	Rated Pressure	0.15 Mpa
	Working Pressure	<0.1Mpa
	Ports	1 x fire ring, 1 x pH, 1 x DO, 1 x PT100, 1 x foam sensor, 1 x exhaust with condenser, spares
	Sterilization	Off-site sterilization (autoclavable)
	Sparge	SS316L ring sparger in the bottom
	Dimension	300×315×708mm
	Net Weight	35 Kg
Agitation	Seal	Top mechanical sealing
	Speed	50 - 1000 rpm
	Blades	3 detachable layers (top: foam breaker; middle: marine blade impeller; bottom: marine blade impeller)
	Baffles	4pcs
Upgrades	Vessel Type	· Single-layer vessel with electric blanket heating; · Jacketed vessel with water (jacket) heating
	Drive	Top Magnetic drive
	Blades	Customizable blades from different options (Spin filter, Cell-lift impeller, Basket impeller, Pitched Blade Impeller, Marine Blade Impeller)

Trust us with your projects!

Lab1st has produced and sold more than 11,000 bioreactors, including 4,500+ Benchtop bioreactors and 6,000+ stainless steel fermenters since 1998. [3 production centers, 20+ years experience]